

**RGDS Stock**  
(Greg A. Perry, Ph.D.)

*RGDS is an inhibitor for PAC-1 binding to activated platelets and is used in platelet activation assays as a specificity control.*

**Equipment:**

1 ml Eppendorf tube

**Reagents:**

2 mg	RGDS (Sigma #A9041)
0.2 ml	Sterile PBS

**Method:**

- 1) Reconstitute the RGDS with the sterile PBS in an Eppendorf tube.

Final stock concentration of RGDS = 10 mg/ml